

Esco Swift™ Extract Automated Nucleic Acid Extraction System



PRINCIPLE

Nucleic acids such as DNA or RNA bind to microscopic magnetic beads found in the accompanying nucleic extraction kit*. A series of steps is involved to (1) lyse the cells containing the nucleic acids, (2) wash away impurities, and (3) release the extracted DNA or RNA from being bound to the magnetic beads in the kit.

**for use with compatible extraction kit*

KEY FEATURES

High Extraction Efficiency

Collection efficiency of up to 98% compared to average market efficiency of 95%.

High Temperature Accuracy

Deep-well heat block accurate heating of wells for effective lysing of cells.

Safe for use

8-strip consumable for placement on rods to prevent cross-contamination and UV light-enabled for decontamination.

User-friendly UI

Real-time display of running program for easy monitoring of extraction protocol.

Versatile blocks and rods

Blocks allow for upgrading to higher throughput.

Standard programs

Capable of storing programs for easy setup of routing protocols.

SPECIFICATIONS

Swift™ Extract, Automated Nucleic Acid Extraction System	
Sample Capacity	32 samples
Sample volume	20-1000 µl
Applicable Consumables	96 deep well plate + 8-tube strip
Uniformity Yield of Wells	CV < 3%
Extraction efficiency	≥ 98%
Operating Temperature	10-40°C
Heating Temperature Control	Lysis and Elution step: 5-125°C
Mixing step	Multiple modes, multiple speed, adjustable
Disinfection / Decontamination	UV Light
Reagent Type	Magnetic bead-based
Communication Interface	USB compatible
Power Supply	100-240 V, 50/60 Hz, 600 W
Dimension (L x W x H)	430 x 395 x 435 mm
Weight	32.5 kg

ORDERING INFORMATION

PRODUCT	Swift™ Extract
Item Code	2210040
Model	SWT-EXT-32, 100-240 VAC
Description	Swift™ Extract, Automated Nucleic Acid Extraction

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Саранск (8342)22-96-24
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

eus@nt-rt.ru || <https://escolifesciences.nt-rt.ru>